S02-500X Digital Microscope User Manual

([English](#))

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Thank you for Purchasing our S02-500X Digital Microscope, it is a slim type tool and can be easily used in different fields including

1. Textile industrial for textile inspection
2. Printing inspection
3. Industrial inspection: PCB, Precision machinery...
4. Educational purpose
5. Hair examination
6. Skin examination
7. Microbiological observation
8. Jewelry & coin (Collections) inspection
9. Visual Assistance
10. Others

Currently we have two models of Digital Microscope: USB port and TV port formats.

The USB microscope must be connected to the computer and TV microscope be connected to visual device (N/P). User interface is very friendly and you can do the Measure work easily.

You can find more information of the product on this manual.

CONTENTS

Introduction	30
Parts of Microscope	31
Specification	32
Notes before use	33
Product Outlook & Standard parts	33
How to install the Driver software	35
How to install the Hotviewer software	38
1X-2X detection instructions	51

Parts of Microscope

- ① LED Light ② Lens ③ Magnification 1X~500X ④ Focus roller
⑤ SNAP & Zoom switch ⑥ Body ⑦ USB line

Specification

Image CMOS Sensor

Controller High Speed DSP (Driver Free available)

Focus Range 0mm ~ 40mm

Snap Shot Software and Hardware

Video Capture Resolution 0.3M.

Built-in 8 White-light LED and adjustable illumination ensure the magnified images are clear and bright

Still Image Capture Resolution 320*240, 640*480, 800*600,
1280*960,1600*1200

Frame Rate 30 f/s under 600 LUX Brightness

Digital Zoom 5X Sequence Mode

Brightness Control Manual adjustment

Magnification Range 1X ~ 500X (Manually)

Power Supply USB Port (5V DC)

USB 2.0 & USB 1.1 Compatible

Operation System Windows XP, Vista, Win 7,Win 8,Win 10

Language: English, Chinese and other language can be OEM

System Requirement: Pentium Computer with 700M Hz & above,
20M HD Space CD ROM Driver, 64MB RAM, Direct X VGA Card

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Product net weight: 320g

Available color: Black, and OEM other colors.

Notes before use

1. Don't disassemble the digital Microscope or change the interior parts, it can cause damage.
2. Don't clean the Microscope with alcohol organic solvents
3. Don't touch the lens with your fingers.
4. Avoid outdoor use if possible.
5. Storage temperature , 0°C ~ 40°C, Humidity: 45%RH ~ 85RH%.
6. In case the product gets wet, leave PC connection immediately. And do not disassemble or dry by hair dryer. Send to repair center if the digit microscope was effected by liquid or other elements.
7. Measurement Data only for reference.

Product Outlook & Standard parts

1. Digital Microscope (1pcs)
2. Calibration ruler (1pcs)
3. Lift Stand (1pcs)

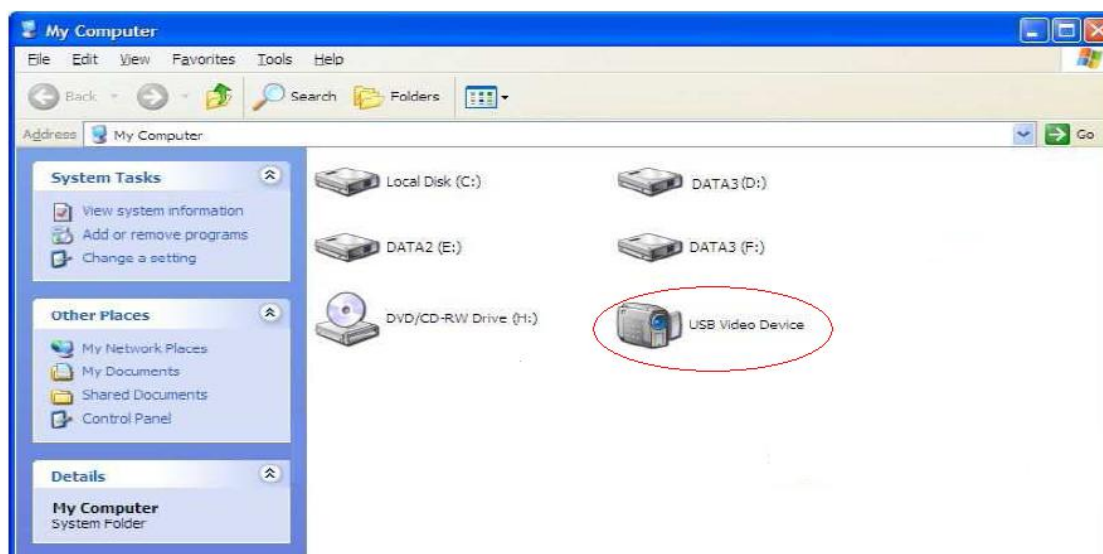
Hardware system requirements

Windows 2000、XP、VISTA、WIN7 Pentium 1G, Celeron, AMD 1G & above, 128MB Memory, 150MB Hard Disc memory space, 16-bit & above VGA ,CD-ROM, USB2.0 or USB1.1.

How to install the Driver

S02-500X Digital Microscope is a free Driver Product, it connected to Windows XP system, it can work immediately.

1. Connect the Portable digital microscope by USB2.0 cable to your computer, double click 'my computer' icon on your computer desktop, below interface will be shown up.

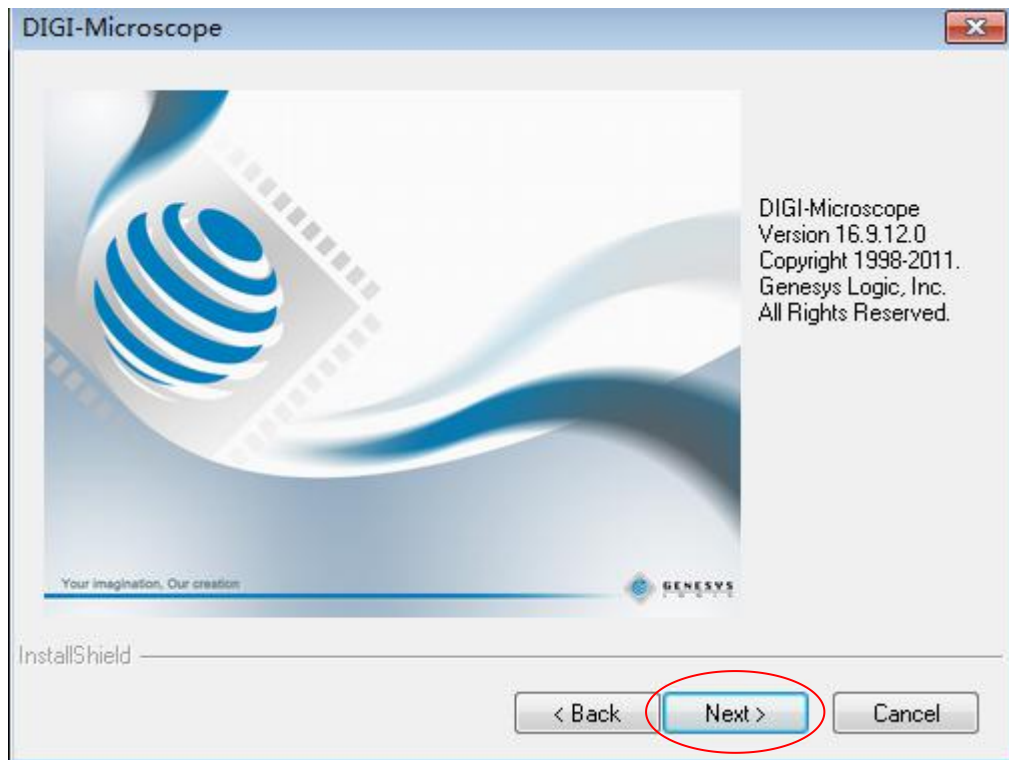
2. Double click the **Video device icon** as above red marked in your computer. Now it is ready to use.

1. Driver installation

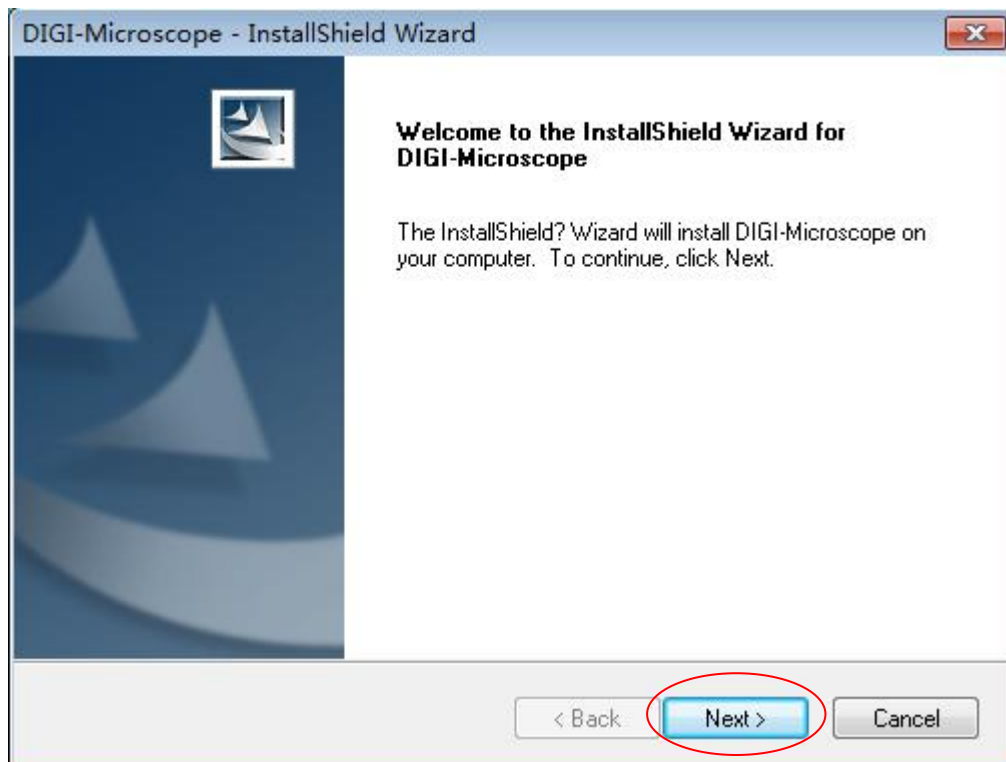
In the computer browser site bar input link: www.indmu.com/hot/2.rar, download the file after extracting, and then double-click the installation software Driver Steup.exe

1 **Connect the product to your computer first**, then double click


 Drive Setup Driver Steup.exe to continue, the screen will show the installation progress.


Click "next" to continue


Click“next”to continue

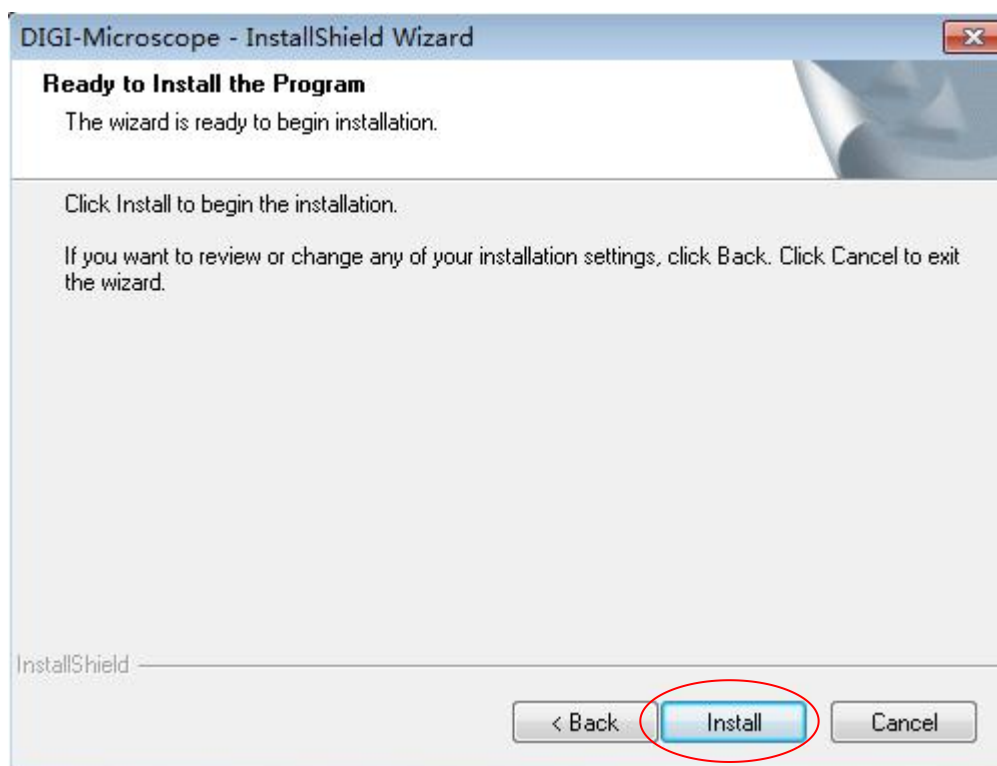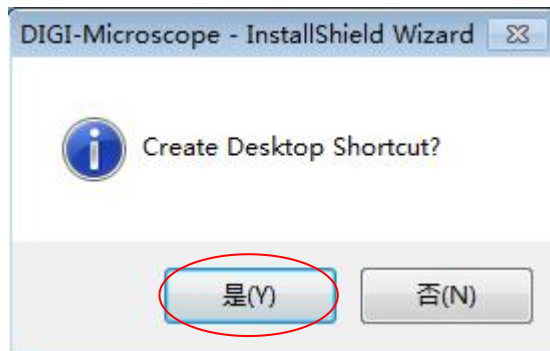Click“Install”to continue, as followings:


Click“是”to continue, as followings:

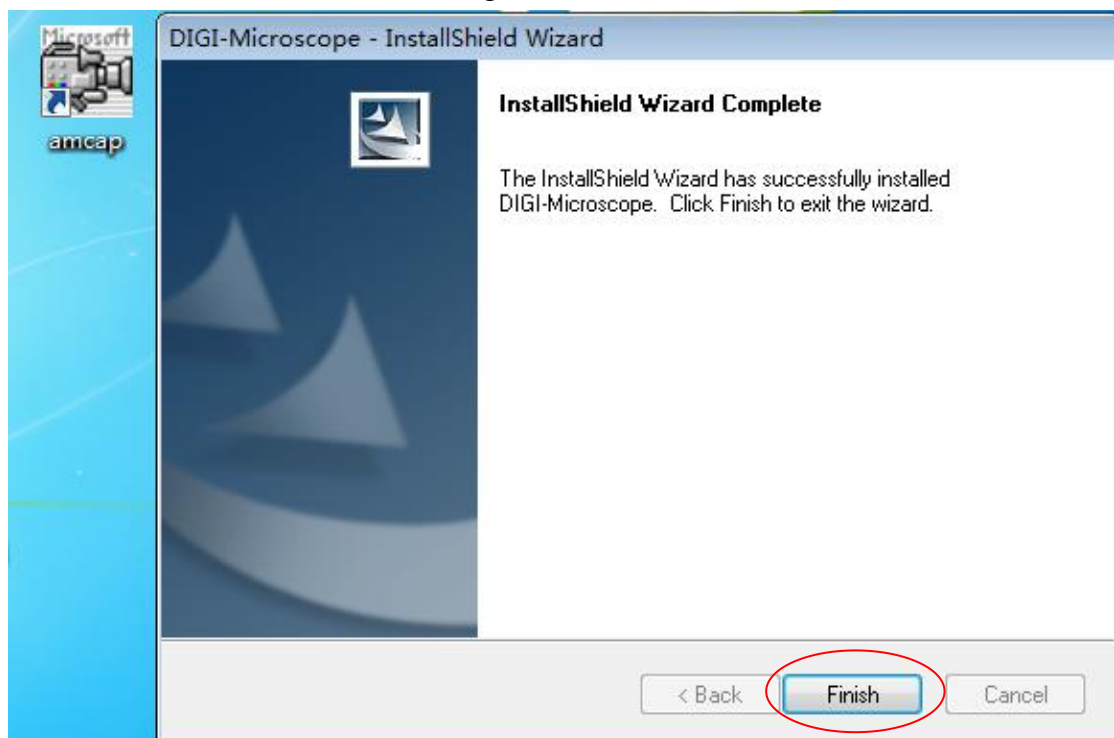Click“Finish”to continue, as followings

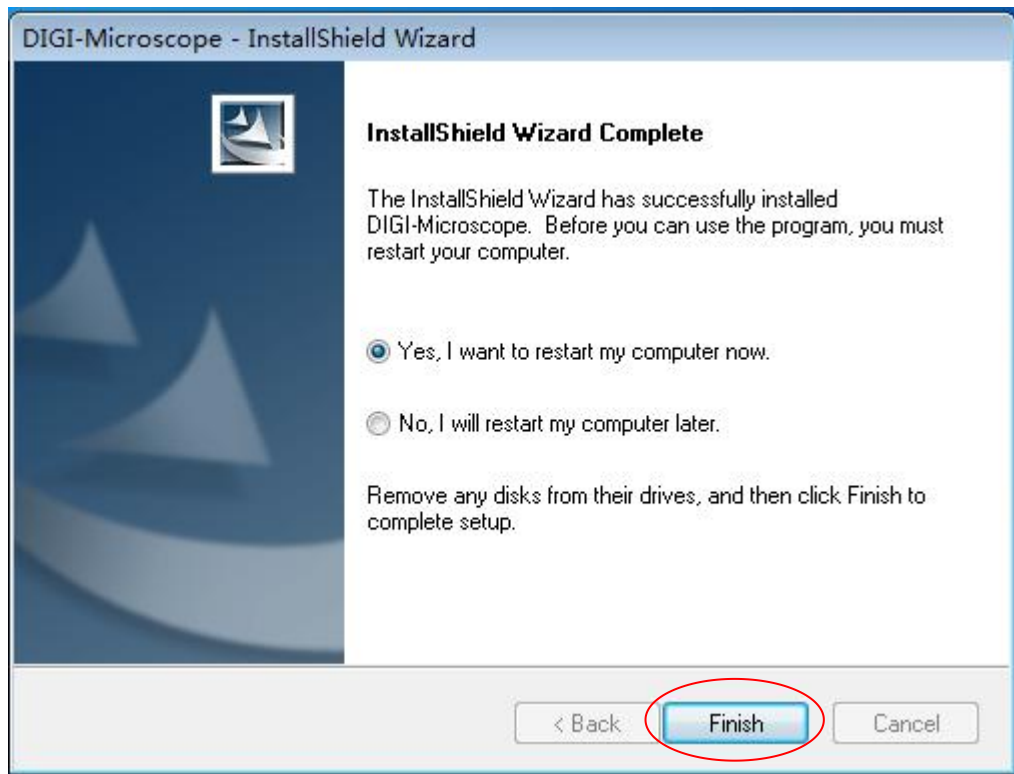Select 'Yes, I want to restart my computer now', and click “finish”.

After restart your computer, Plug-in the USB port of Digital Microscope into Computer USB port. Following the user guide by steps and complete by click “Next” -> “Continue anyway” -> “Finish”.

2 Hotviewer installations: **Windows(PC)** (Win7 for example)

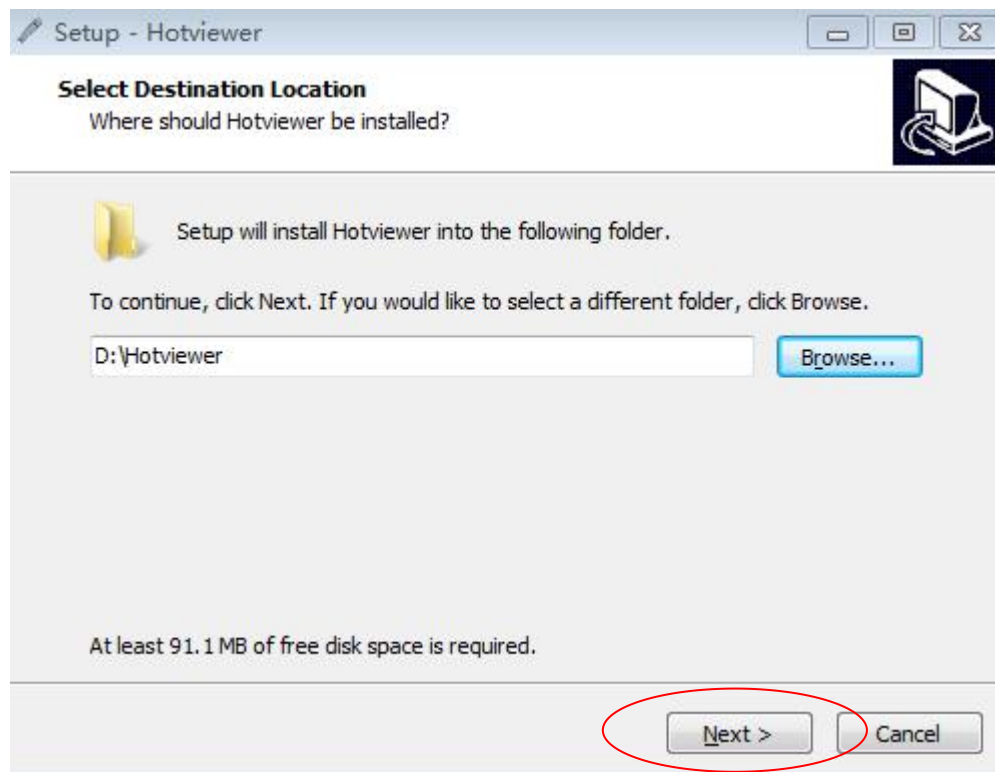
Reminder: When installing Hotviewer software, please exit the anti-virus software to install, because anti-virus software will hinder the installation

process and destruction of software, installation errors and can cause an unsuccessful installation and other anomalies.


1. In the computer browser site bar input link: www.indmu.com/hot/2.rar, download the file after extracting, and then double-click the installation software Hotviewer.exe


Click "OK"


Click "Next"

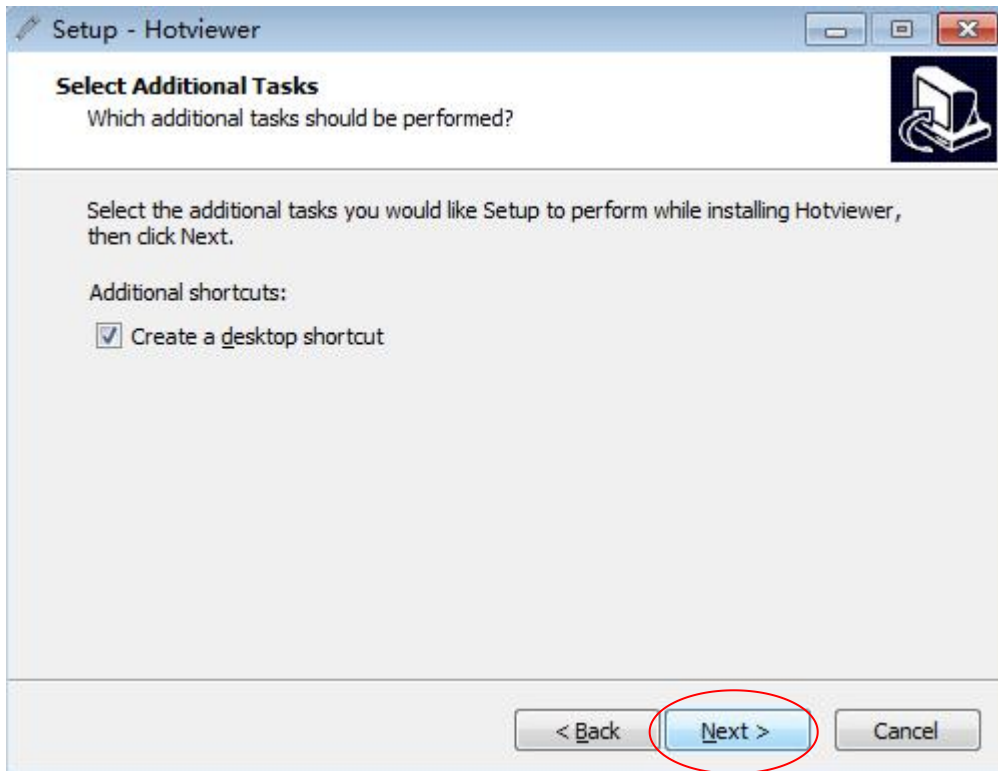
Click “Next”

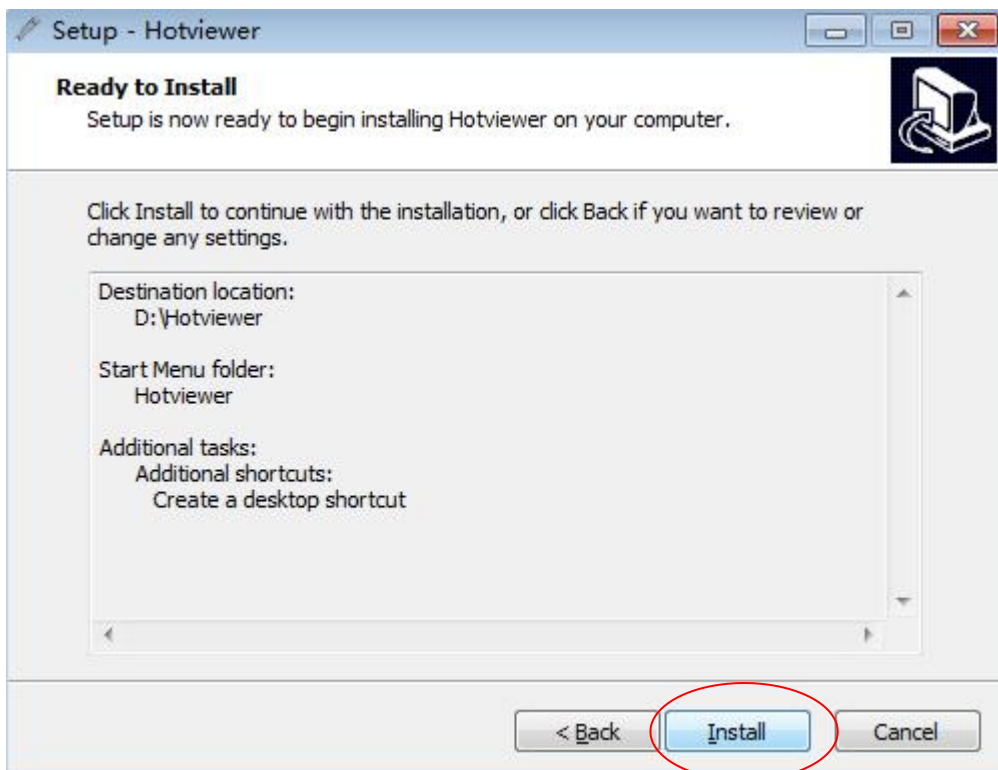Click “Install”

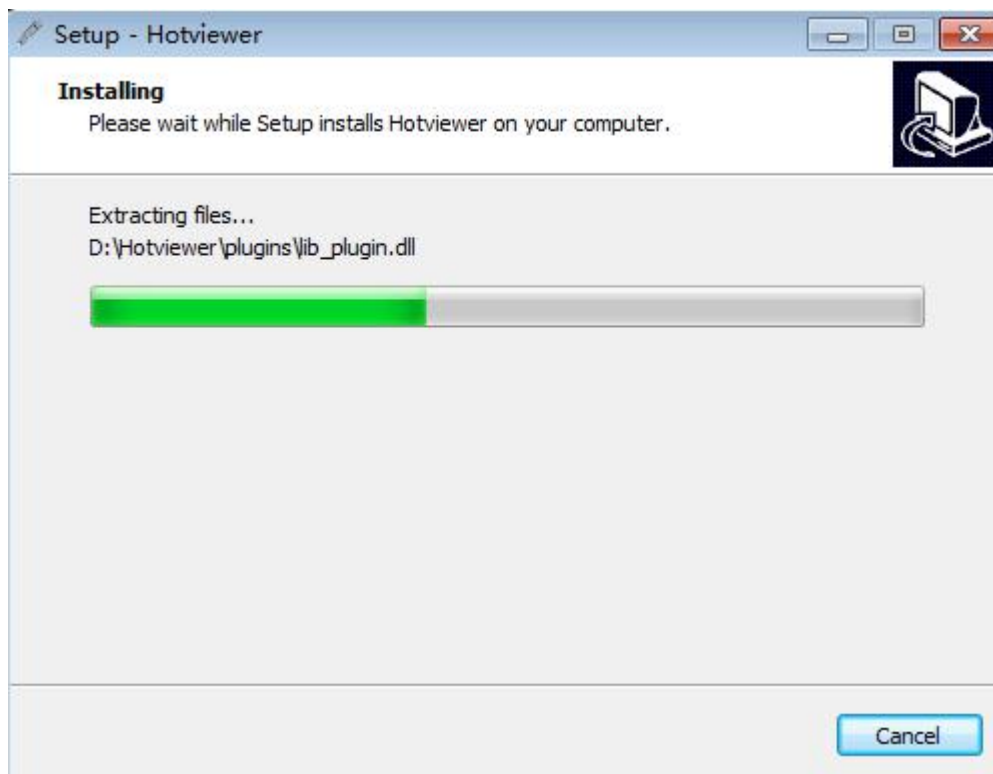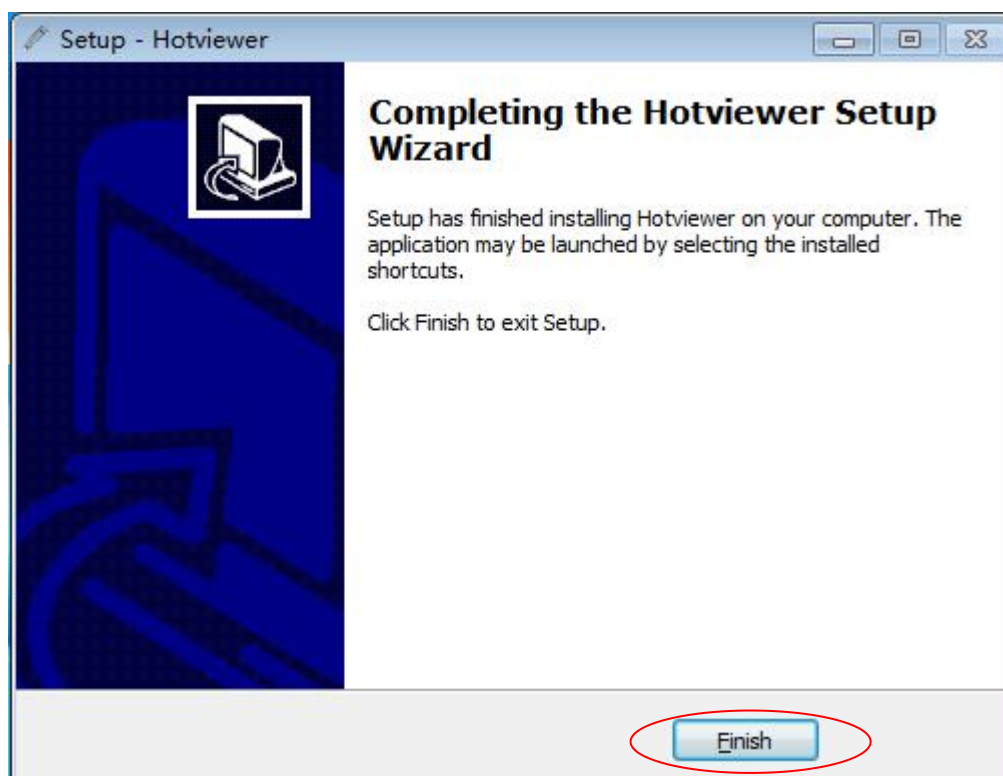Waiting for the installation progress to complete


Click "Finish" to complete the software installation

2. When a microscope was connected to a computer via a USB cable,

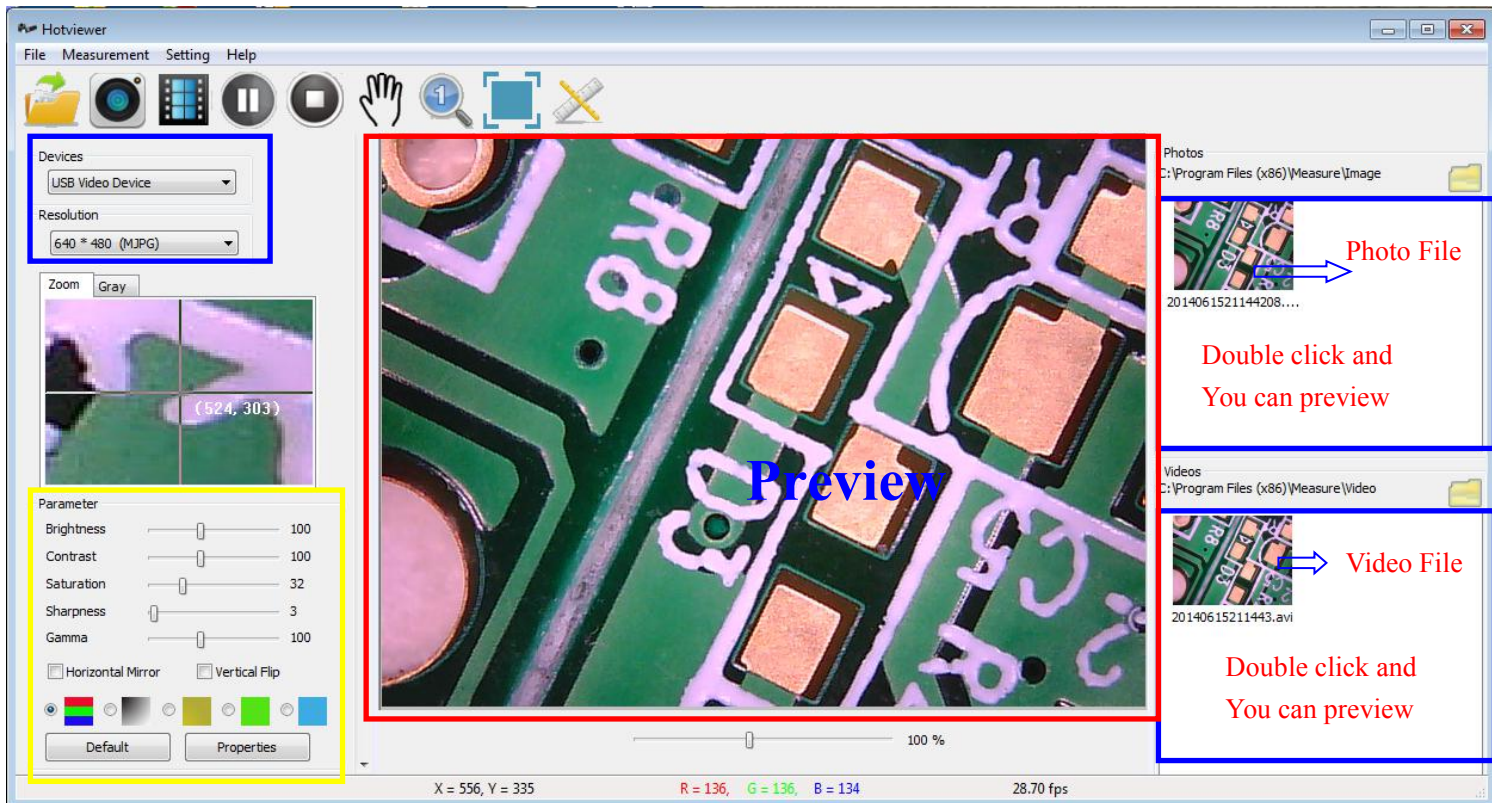Click the **Hotviewer**, you can open the preview of the application interface.

3. Interface introduction

1. Tool bar

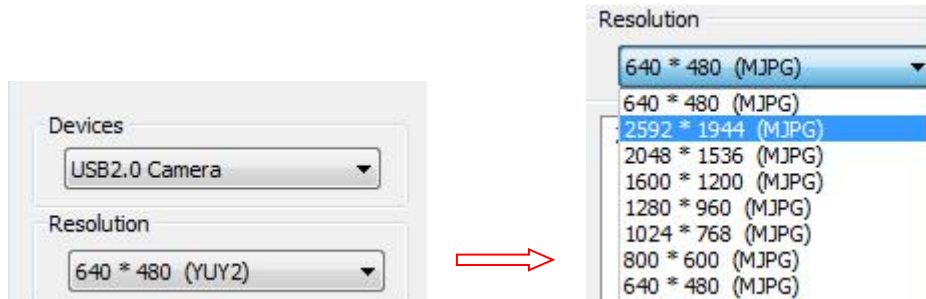
Open image **photo** video **pause/play** stop


move initial full screen measurement


2. Devices

Digital zoom

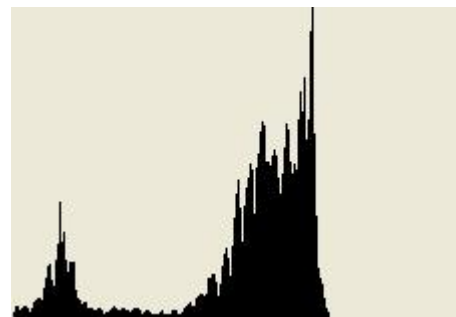Gray level

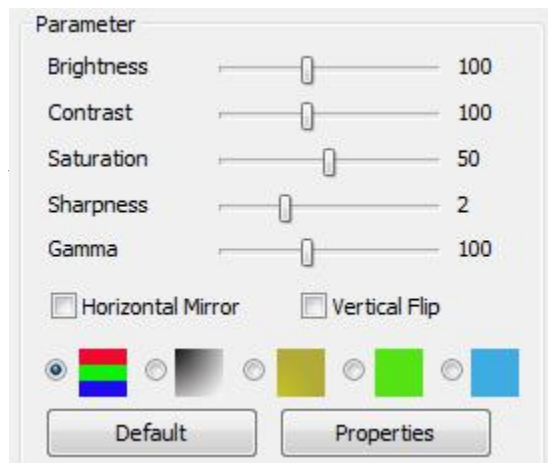
You can set the parameter, if you want to restore the factory state, please Click 'Default'.


3.menu

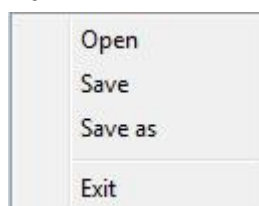File


Measurement


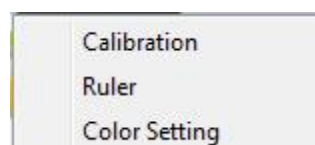
Click this menu and open the measurement interface.

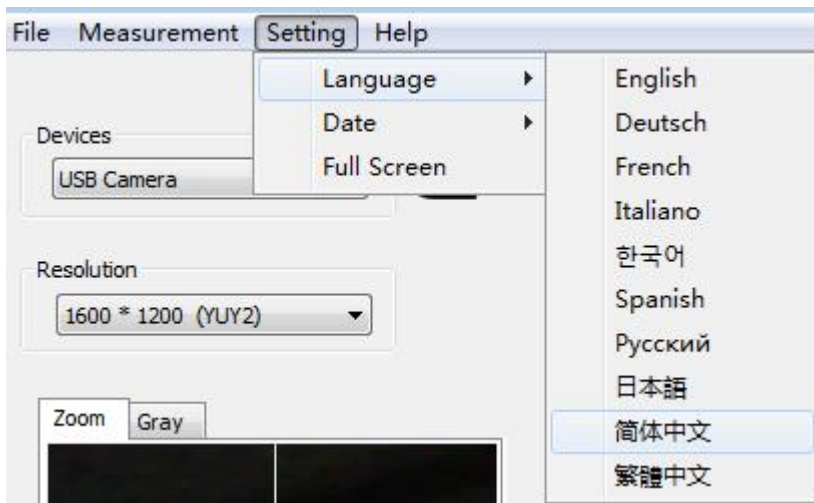Setting

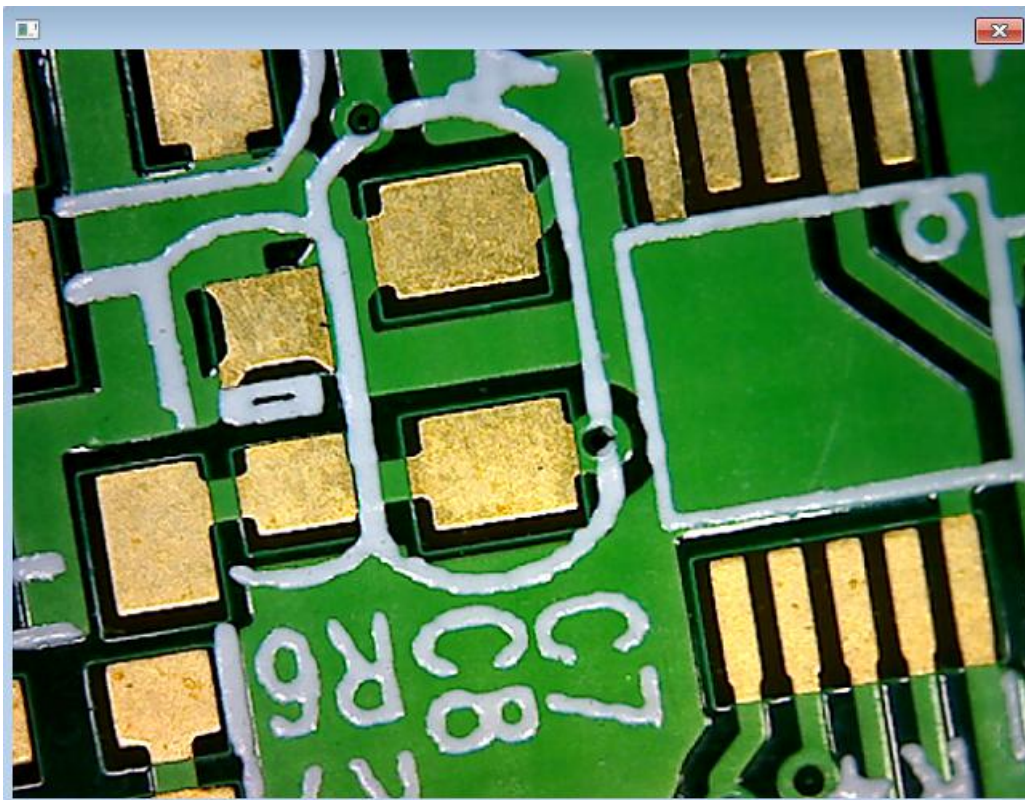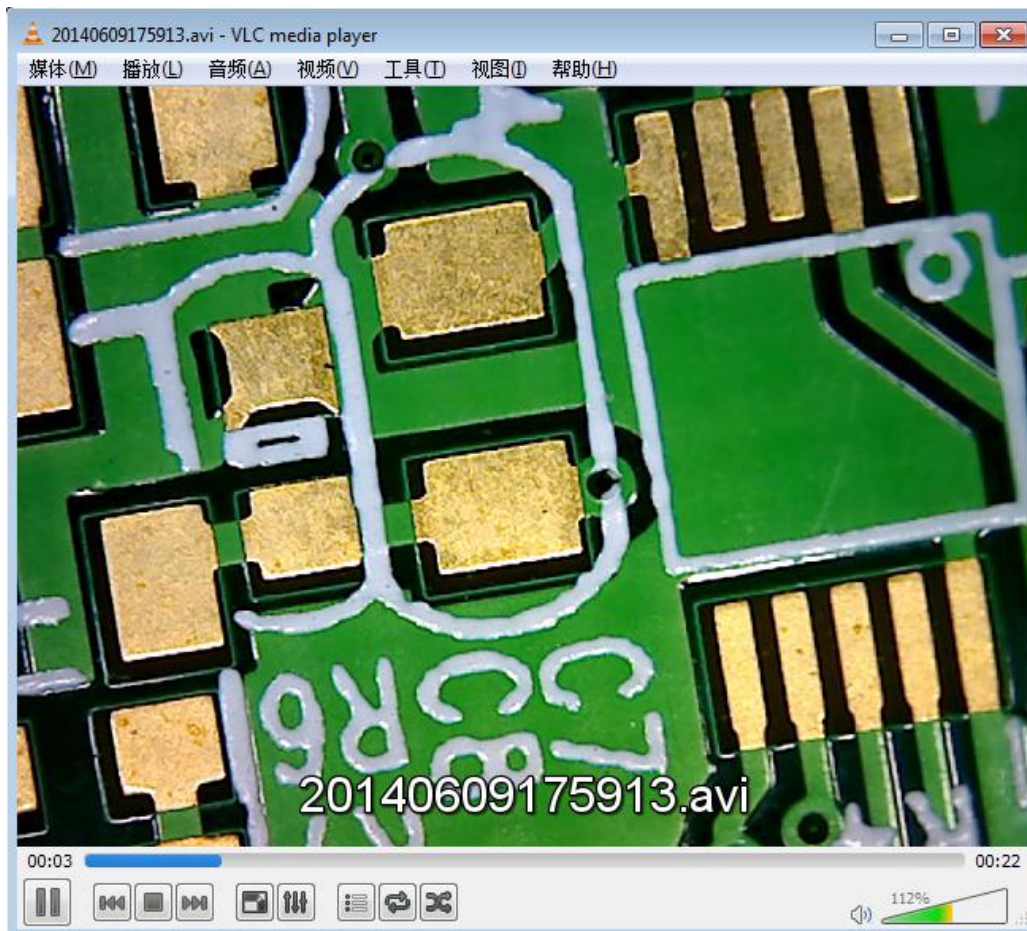
4 .Photo & Video (double click file and open it)

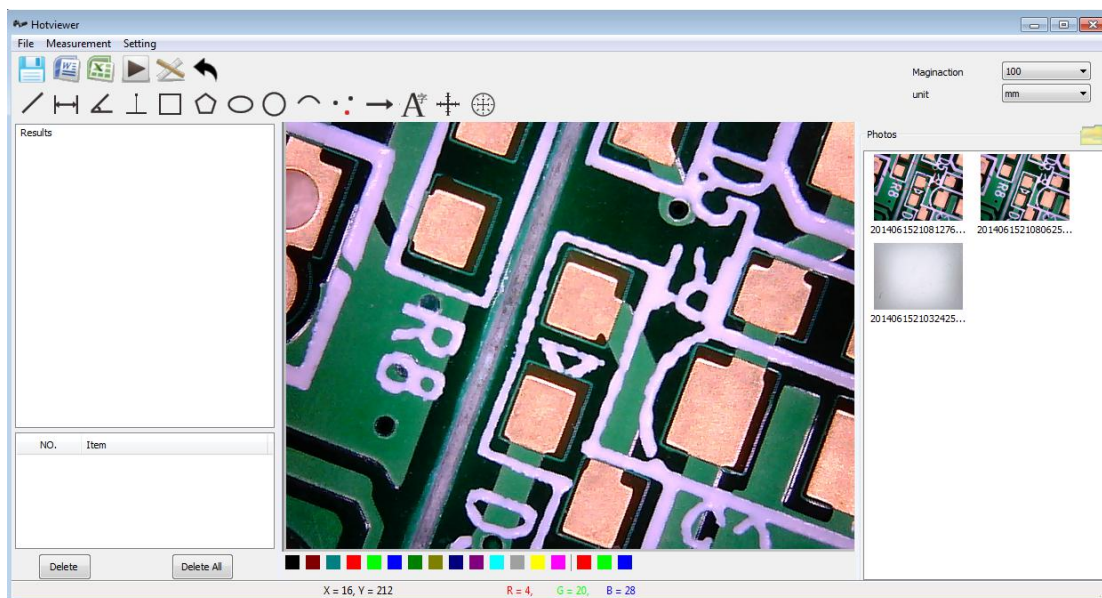
5. Measurement interface

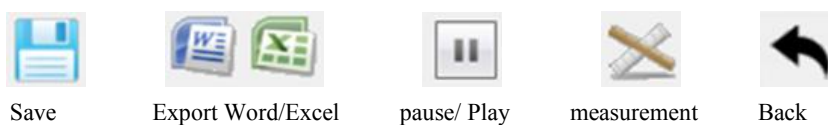6. Tool bar


Save

Export Word/Excel

pause/ Play


measurement

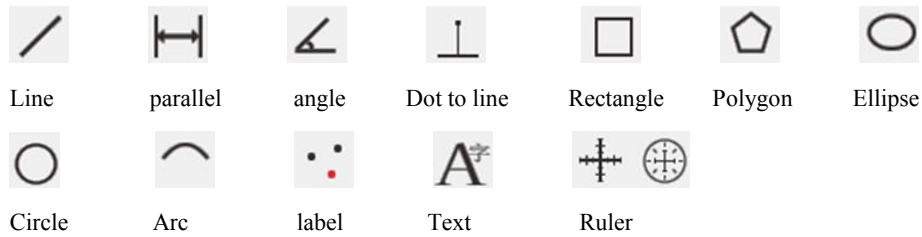Back


7. menu introduction

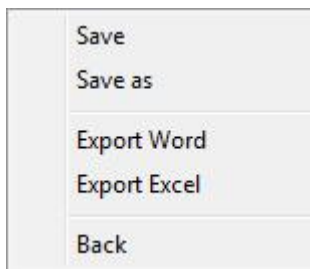File

Measurement

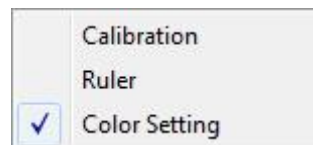Setting

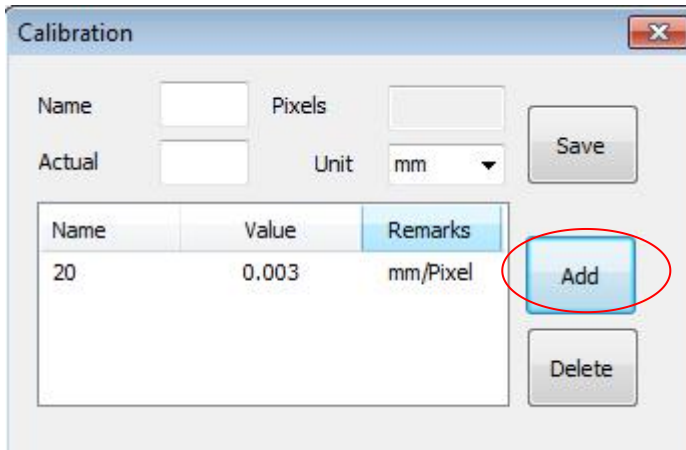
8. Measurement


(1) Calibration

Before measurement, please do calibration to make sure a certain accuracy. Our experience tells us that the measurement accuracy can be controlled to as high as 99%. Human eye distortion or defocus could cause measurement tolerance. More practice is helpful to measurements.


Step 1 ,Click , and press ' add'

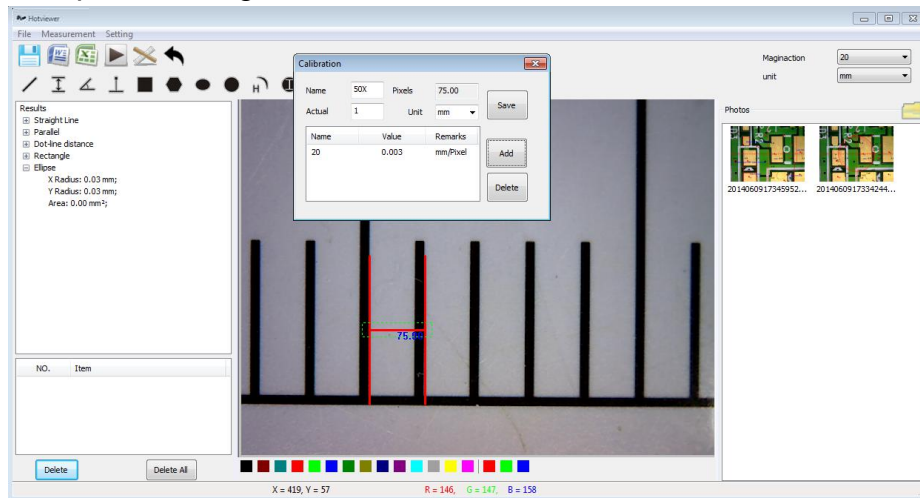
Step2, Shooting calibration ruler calibrate the ruler

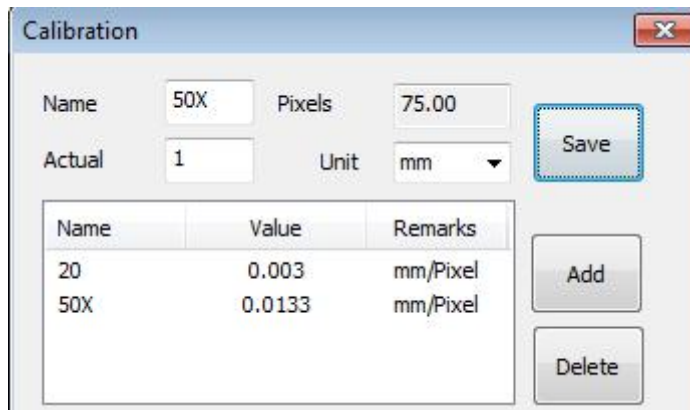
Step3, Record key information and save it

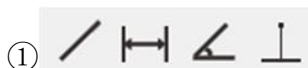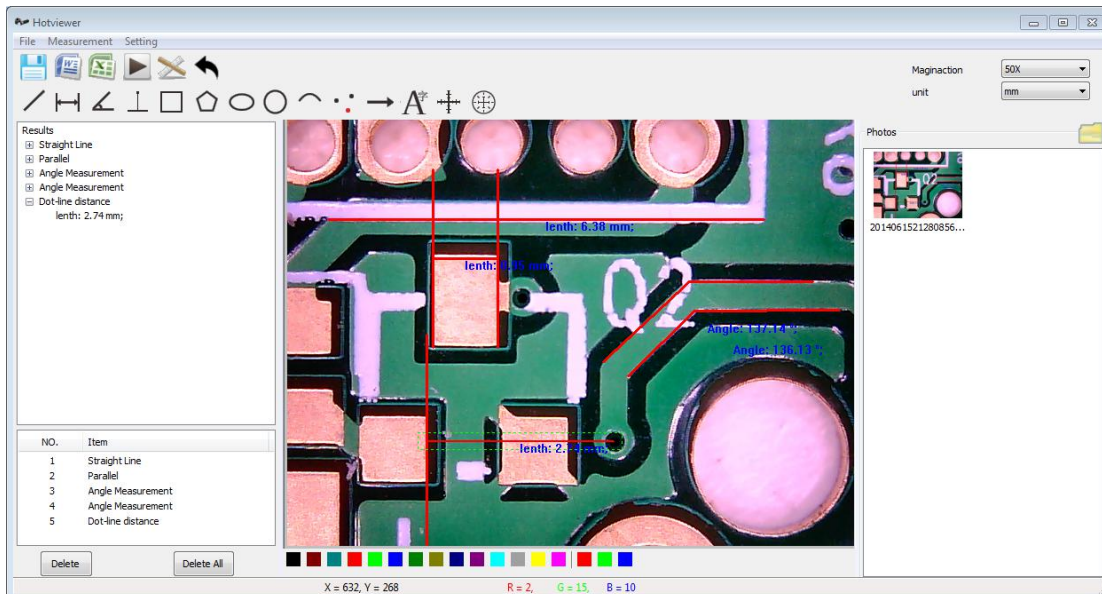
2. Measurement


Press  to save picture.

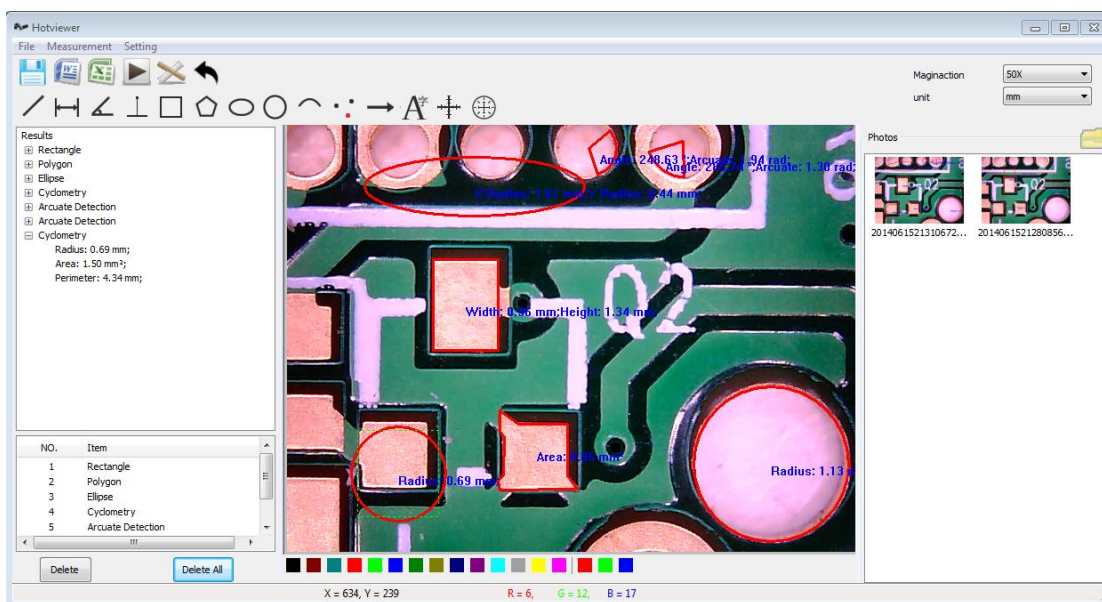Press  to save picture.

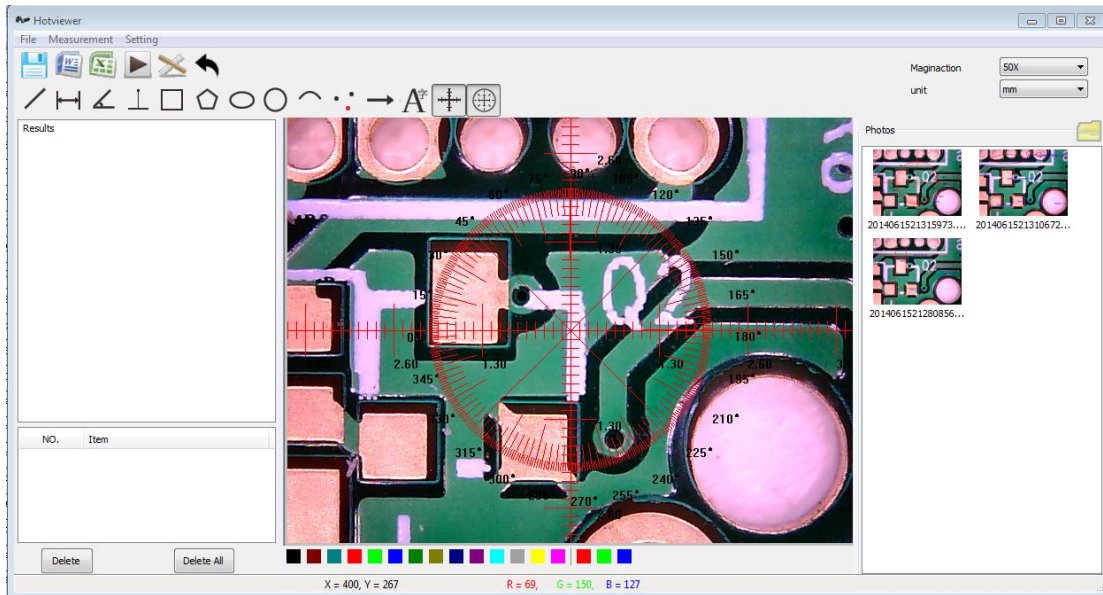Press  to save picture.

④ Export result

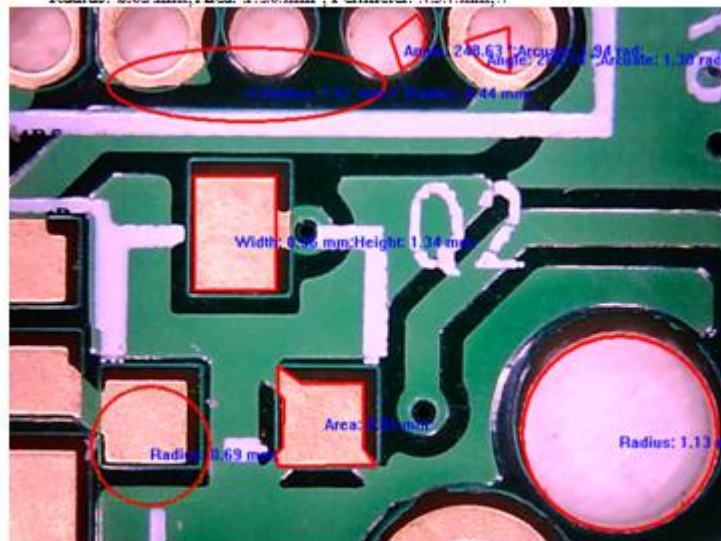
Press , then export the file selected.

Item: 20140615213106723.bmp


Data: 2014-06-15 21:33

Magnification ruler 0.0130 mm/px

1. Rectangle:.
Width: 0.96 mm; Height: 1.34 mm; Area: 1.29 mm²;
2. Polygon:.
Area: 0.95 mm²;
3. Ellipse:.
X Radius: 1.61 mm; Y Radius: 0.44 mm; Area: 2.21 mm²;
4. Cyclometry:.
Radius: 1.13 mm; Area: 4.03 mm²; Perimeter: 7.11 mm;
5. Arcuate Detection:.
Angle: 285.74 °; Arcuate: 1.30 rad; Radian: 0.69 mm;
6. Arcuate Detection:.
Angle: 248.63 °; Arcuate: 1.94 rad; Radian: 0.79 mm;
7. Cyclometry:.
Radius: 0.69 mm; Area: 1.50 mm²; Perimeter: 4.34 mm;

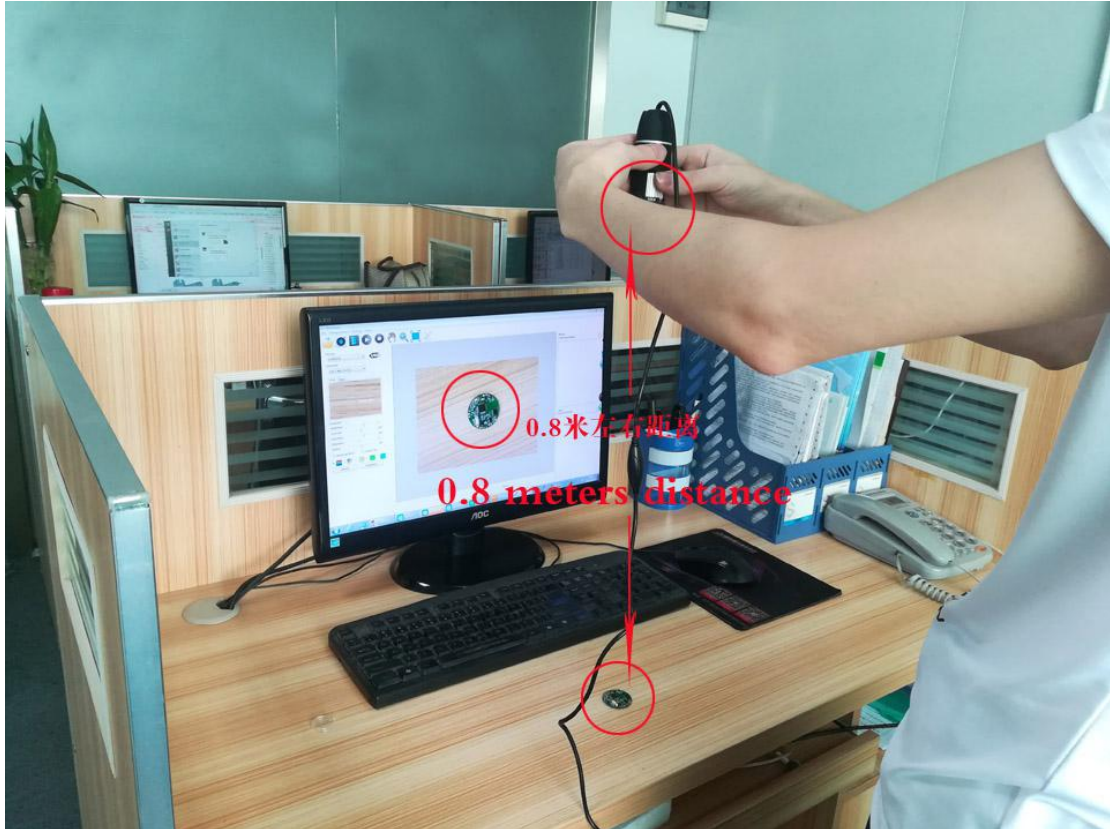
Instructions for detecting objects and zooming 1-2X: when microscope lens at the distance of 0.8 meters or so away from the object, to zoom 1.5-2X, as shown below

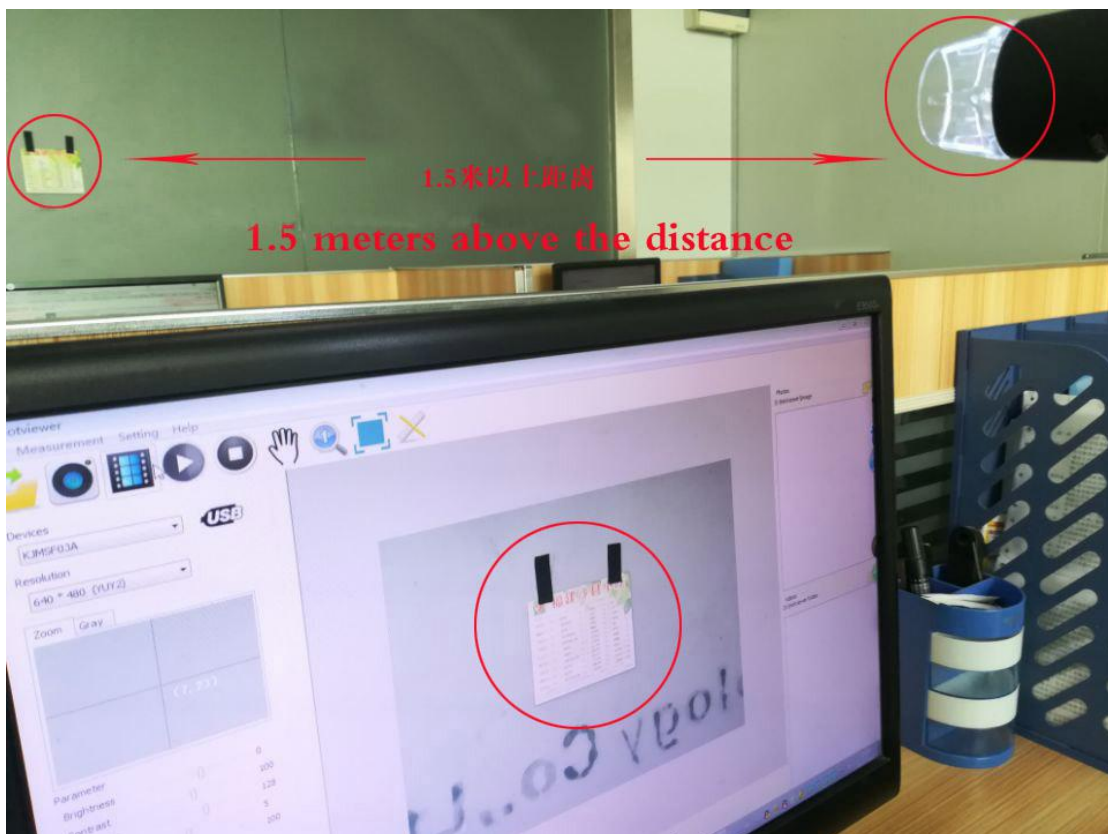when microscope lens at the distance of more than 1.5meters
or so away from the object, to zoom 1X, as shown below

To zoom 1-2X, your hands need to hold the microscope and turn to the 50X scale, fine-tuning focus and detect, when to get a clear image, keep your hands stable. Do not move, if you move the arm slightly, the image will move too, resulting in the image not clear. If you want to make the image more clearly to zoom 1-2 time, you need to keep the microscope from the object more than 0.8 meters away, and fix it unmovable, and then zoom 1-2X detection. On the standard black plastic lift stand, it can not be zoomed 1X.